

July 4, 2026

Laboratory 12

Delete Antibody Data by Project

Excel

19th IHIWS database system development team

Hide or Delete?

Decide before you act. Deleting cannot be undone.

Just hide it? Use Data Visibility

If you only want to stop one project from seeing the data, do NOT delete it. Open the Data Visibility tab and hide it. This is reversible - you can show it again at any time.

Really remove it? Use Delete

Use Delete only to permanently remove the data. It is erased from your lab and disappears from ALL linked projects, not only the one you selected. This cannot be undone.

Delete Antibody Data by Project

<https://data.ihiw19.org>

IHIW19

[User Guide](#)

Dashboard

[Lab View](#)

[Project View](#)

 Doc User ▾

Dashboard

Click "Lab View"

Welcome to IHIW19 Management System, Doc User

Metadata Completion Status

HLA Metadata

Sample Metadata (Stage 1)

Sample: 0 of 0 completed (0%)

Subject Metadata (Stage 2)

Subject: 0 of 0 completed (0%)

No HLA metadata found

Antibody Metadata

Sample Metadata (Stage 1)

Sample: 0 of 3 completed (0%)

Subject Metadata (Stage 2)

Subject: 0 of 3 completed (0%)

No Antibody metadata found

Lab View

Manage lab members, upload files, and configure sharing settings with projects.

[Go to Lab View](#)

Project View

Manage projects, view shared files, and import/export data.

[Go to Project View](#)

Delete Antibody Data by Project

<https://data.ihw19.org/lab>

IHIW19

[User Guide](#)

[Dashboard](#)

Lab View

[Project View](#)

 Doc User ▾

Lab View

Lab: DOCLAB03

Samples not visible to any project: HLA: 0 / Antibody: 0

Samples visible per project:

- HLA Typing Workshop 2025: HLA 0 (hidden: 0) / Antibody 0 (hidden: 0)
- Antibody Screening 2025: HLA 0 (hidden: 0) / Antibody 0 (hidden: 0)

 Lab Members

 Project Linkage

 Genotype Data

 Antibody Testing Data

 Sample Metadata

 Subject Metadata

 Data Visibility

 Genotype Sample Visibility

 Token Management

Click “Antibody Testing Data”

Lab Members

 Add Lab Member

Name	Email	Role	Status	Action
Doc User	pi@doclab03.example.com	PI	Active	

Delete Antibody Data by Project

<https://data.ihw19.org/lab/#antibody-testing-data>

IHIW19

User Guide

Dashboard

Lab View

Project View

Doc User

Lab View

Lab: DOCLAB03

Samples not visible to any project: HLA: 0 / Antibody: 0

Samples visible per project:

- HLA Typing Workshop 2025: HLA 0 (hidden: 0) / Antibody 0 (hidden: 0)
- Antibody Screening 2025: HLA 0 (hidden: 0) / Antibody 0 (hidden: 0)

Antibody Testing Data list

Lab Members

Project Linkage

Genotype Data

Antibody Testing Data

Sample Metadata

Subject Metadata

Data Visibility

Genotype Sample Visibility

Token Management

Antibody Testing Data Management

Import Antibody Data

Delete Selected

Guide

Project

Visibility

Search

Items per page

All linked projects

All

Search...

10



Sample ID

Specificity

Visibility

Actions



SAMP_001

A*01:01, A*02:01

—



DELETE



SAMP_002

B*07:02

—



DELETE



SAMP_003

DRB1*15:01

—



DELETE

Delete Antibody Data by Project

<https://data.ihw19.org/lab/#antibody-testing-data>

IHIW19

User Guide

Dashboard

Lab View

Project View

Doc User

Lab View

Lab: DOCLAB03

Select a linked project. The list now shows only what that project can see (Visibility = Visible)

Samples not visible to any project: HLA: 0 / Antibody: 0

Samples visible per project:

- HLA Typing Workshop 2025: HLA 0 (hidden: 0) / Antibody 0 (h
- Antibody Screening 2025: HLA 0 (hidden: 0) / Antibody 0 (hid

Lab Members

Project Linkage

Genotype Data

Antibody Testing Data

Sample Metadata

Subject Metadata

Data Visibility

Genotype Sample Visibility

Token Management

Antibody Testing Data Management

Import Antibody Data

Delete Selected

Guide

Project

Visibility

Search

Items per page

HLA Typing Works

Visible

Search...

10

Filter active: showing samples VISIBLE to "HLA Typing Workshop 2025" (2 shown). Owner view is filtered.

Clear filter

☐	Sample ID	Specificity	Visibility	Actions
--------------------------	-----------	-------------	------------	---------

☐	SAMP_001	A*01:01, A*02:01	Visible	 DELETE
--------------------------	----------	------------------	---------	--

☐	SAMP_003	DRB1*15:01	Visible	 DELETE
--------------------------	----------	------------	---------	--

Delete Antibody Data by Project

<https://data.ihw19.org/lab/#antibody-testing-data>

IHIW19

User Guide

Dashboard

Lab View

Project View

Doc User

Lab View

Lab: DOCLAB03

Samples not visible to any project: HLA: 0 / Antibody: 0

Samples visible per project:

- HLA Typing Workshop 2025: HLA 0 (hidden: 0) / Antibody 0 (hidden: 0)
- Antibody Screening 2025: HLA 0 (hidden: 0) / Antibody 0 (hidden: 0)

Visibility filter: All / Visible / Hidden

Each row shows a badge:
Visible = the project can see it
Hidden = hidden from the project

Lab Members

Project Linkage

Genotype Data

Antibody Testing Data

Sample Metadata

Subject Metadata

Data Visibility

Genotype Sample Visibility

Antibody Testing Data Management

Antibody Data

Delete Selected

Guide

Project

Visibility

Search

Items per page

HLA Typing Works

All

Search...

10



Sample ID

Specificity

Visibility

Actions



SAMP_001

A*01:01, A*02:01

Visible



DELETE



SAMP_002

B*07:02

Hidden



DELETE



SAMP_003

DRB1*15:01

Visible



DELETE

Choose "All" to see everything you own (Visible and Hidden): the badge tells you how each

Delete Antibody Data by Project

<https://data.ihw19.org/lab/#antibody-testing-data>

IHIW19

[User Guide](#)

[Dashboard](#)

[Lab View](#)

[Project View](#)

[Doc User](#) ▾

Lab View

Lab: DOCLAB03

A banner shows the active filter. Click "Clear filter" to see everything again

Set Visibility = Hidden to review only the samples hidden from this project

Samples not visible to any project: HLA: 0 / Antibody: 0

Samples visible per project:

- HLA Typing Workshop 2025: HLA 0 (hidden: 0) / Antibody 0 (hidden: 0)
- Antibody Screening 2025: HLA 0 (hidden: 0) / Antibody 0 (hidden: 0)

 Lab Members

 Project Linkage

 Genotype Data

 **Antibody Testing Data**

 Sample Metadata

 Subject Metadata

 Data Visibility

 Genotype Sample Visibility

 Token Management

Antibody Testing Data Management

[Import Antibody Data](#)

[Delete Selected](#)

[Guide](#)

Project

Visibility

Search

Items per page

HLA Typing Works ▾

Hidden ▾

Search...

10

Filter active: showing samples HIDDEN from "HLA Typing Workshop 2025" (1 shown). Owner view is filtered.

[Clear filter](#)



Sample ID

Specificity

Visibility

Actions



SAMP_002

B*07:02

Hidden



DELETE

Delete Antibody Data by Project

<https://data.ihw19.org/lab/#antibody-testing-data>

IHIW19 User Guide Dashboard **Lab View** Project View

Doc User ▾

Lab View

Lab: DOCLAB03

Samples not visible to any project: HLA: 0 / Antibody: 0

Samples visible per project:

- HLA Typing Workshop 2025: HLA 0 (hidden: 0) / Antibody 0 (hidden: 0)
- Antibody Screening 2025: HLA 0 (hidden: 0) / Antibody 0 (hidden: 0)

Samples filtered out (not shown) are NOT deleted

Lab Members

Antibody Testing Data Management

Import Antibody Data

Delete Selected

Guide

Delete Selected removes only the rows shown here and checked

Antibody Testing Data

Sample Metadata

Subject Metadata

Data Visibility

Genotype Sample Visibility

Search

Items per page

Search...

10

Filter: showing samples HIDDEN from "HLA Typing Workshop 2025" (1 shown). Owner is filtered. Clear filter



Sample ID

Specificity

Visibility

Actions



SAMP_002

B*07:02

Hidden



DELETE

But the delete is still permanent and lab-wide: the shown samples are removed from ALL

Delete Antibody Data by Project

<https://data.ihw19.org/lab/#antibody-testing-data>

IHW19

User Guide

Dashboard

Lab View

Project View

Doc User

Lab View

Lab: DOCLAB03

Samples not visible to any project: HLA: 0 / Antibody: 0

Samples visible per project:

- HLA Typing Workshop 2025: HLA 0 (hidden: 0) / Antibody 0 (hidden: 0)
- Antibody Screening 2025: HLA 0 (hidden: 0) / Antibody 0 (hidden: 0)

Then click "Delete Selected"

Check the rows to delete

Lab Members

Project Linkage

Genotype Data

Antibody Testing Data

Sample Metadata

Subject Metadata

Data Visibility

Genotype Sample Visibility

Token Management

Antibody Testing Data Management

Import Antibody DataDelete SelectedGuide

Project: HLA Typing Works ▼Visibility: All ▼Search: Search... 🔍Items per page: 10 ▼

Sample ID	Specificity	Visibility	Actions
☒ SAMP_001	A*01:01, A*02:01	Visible	☒ ☐ DELETE
☒ SAMP_002	B*07:02	Hidden	☒ ☐ DELETE
☒ SAMP_003	DRB1*15:01	Visible	☒ ☐ DELETE

Delete Antibody Data by Project

IHIW1 Doc User

Lab View

Lab: DOCLAB03

Samples not visible to any project: HLA: 0 / 0

Samples visible per project:

- HLA Typing Workshop 2025: HLA 0 (hidden: 0)
- Antibody Screening 2025: HLA 0 (hidden: 0)

Lab Members

Project Linkage

Antibody Data Delete Selected Guide

Items per page 10

Sample ID	Specificity	Visibility	Actions
SAMP_001	A*01:01, A*02:01	Visible	DELETE
SAMP_002	B*07:02	Hidden	DELETE
SAMP_003	DRB1*15:01	Visible	DELETE

Warning: the delete removes the data from ALL linked projects

Are you sure you want to delete the selected antibody data?

Selected antibody data.

Warning: This permanently removes the data from your lab. It will disappear from **all linked projects**, not only the selected one. This action cannot be undone.

3 of 3 selected sample(s) are also visible to other linked projects.
[Show details](#)

Cancel Delete

Impact summary: how many selected samples other projects still see

Sample Metadata

Subject Metadata

Data Visibility

Genotype Sample Visibility

Token Management

Delete Antibody Data by Project

IHIW19

User Guide

Dashboard

Lab View

Project View

Doc User

Lab View

Lab: DOCLAB03

Samples not visible to any project: HLA: 0 / 0

Samples visible per project:

- HLA Typing Workshop 2025: HLA 0 (hidden: 0)
- Antibody Screening 2025: HLA 0 (hidden: 0)

Lab Members

Project Linkage

Genotype Data

Antibody Testing Data

Sample Metadata

Subject Metadata

Token Management

Confirm Deletion



Are you sure you want to delete the selected antibody data?

Selected antibody data: 3

Warning: This permanently removes the data from your lab. It will disappear from **all linked projects**, not only the selected one. This action cannot be undone.

3 of 3 selected sample(s) are also visible to other linked projects.

[Show details](#)

- SAMP_001 → Antibody Screening 2025
- SAMP_002 → Antibody Screening 2025
- SAMP_003 → Antibody Screening 2025

Cancel

Delete

Click "Show details" to list, per sample, which other projects can still see it

Then Click "Delete"

Delete Antibody Data by Project

IHIW19

User Guide

Dashboard

Lab View

Project View

Doc User

Lab View

Lab: DOCLAB03

Deletion Result



Successfully deleted 3 antibody data samples.

Deletion completed

OK

Samples not visible to any project: HLA: 0 / Antibody: 0

Samples visible per project:

- HLA Typing Workshop 2025: HLA 0 (hidden: 0) / Antibody 0 (hidden: 0)
- Antibody Screening 2025: HLA 0 (hidden: 0) / Antibody 0 (hidden: 0)

Lab Members

Project Linkage

Genotype Data

Antibody Testing Data

Sample Metadata

Subject Metadata

Data Visibility

Genotype Sample Visibility

Token Management

Antibody Testing Data Management

Import Antibody Data

Delete Selected

Guide

Project

Visibility

Search

Items per page

HLA Typing Works

All

Search...

10



Sample ID

Specificity

Visibility

Actions



SAMP_001

A*01:01, A*02:01

Visible



DELETE



SAMP_002

B*07:02

Hidden



DELETE



SAMP_003

DRB1*15:01

Visible



DELETE

Delete Antibody Data by Project

<https://data.ihw19.org/lab/#antibody-testing-data>

IHIW19

[User Guide](#)

[Dashboard](#)

[Lab View](#)

[Project View](#)

 Doc User ▾

Lab View

Lab: DOCLAB03

Samples not visible to any project: HLA: 0 / Antibody: 0

Samples visible per project:

- HLA Typing Workshop 2025: HLA 0 (hidden: 0) / Antibody 0 (hidden: 0)
- Antibody Screening 2025: HLA 0 (hidden: 0) / Antibody 0 (hidden: 0)

 Lab Members

 Project Linkage

 Genotype Data

 **Antibody Testing Data**

 Sample Metadata

 Subject Metadata

Antibody Testing Data Management

[Import Antibody Data](#)

 [Delete Selected](#)

 [Guide](#)

Project

Visibility

Search

Items per page

HLA Typing Works ▾

All ▾

Search...



10



Sample ID

Specificity

Visibility

Actions

No antibody data found

The delete is permanent and lab-wide: the data is removed from every linked project, not only the selected one. If a sample has no other data (for example, no Genotype/HLA data), its record is removed as well.